



U.S. Corporate Headquarters
400 Valley Rd.
Warrington, PA 18976
1(800) 523-2575 / (215) 343-6484
1(800)343-3291 fax
info@polysciences.com

Polysciences Europe GmbH
Handelsstrasse 3
D-69214 Eppelheim, Germany
+(49) 6221-765767
+(49) 6221-764620 fax
info@polysciences.de

Polysciences Asia-Pacific, Inc.
2F-1, 207 Tunhwa N. Rd.
Taipei, Taiwan 10595
(886) 2 8712 0600
(886) 2 8712 2677 fax
info@polysciences.tw

TECHNICAL DATA SHEET 521

Page 1 of 1

Aqua-Poly/Mount Coverslipping Medium

INTRODUCTION

Aqua-Poly/Mount is a non-permanent aqueous mounting medium formulated for coverslipping directly from aqueous solutions. It is non-fluorescing and has an antifade component to increase the viewing time of the specimen. Use Aqua-Poly/Mount with most fluorescent dyes and stains including DAB, Alkaline Phosphatase Fast Red, AEC (aminoethylcarbazole) and a variety of other chromogens to enhance and retain fluorescent intensity.

Aqua-Poly/Mount can be used for frozen sections, fat stains, chromogens for immunohistochemistry and in situ hybridization as well as other applications requiring a water soluble mounting medium. Using Aqua-Poly/Mount, the coverslip will dry in 24 to 48 hours depending on the amount of medium used. **Aqua-Poly/Mount solidifies under coverslip on the microscope slide after 24 hours with the following laboratory conditions and ambient temperature; relative humidity in winter time at greater than 20% and relative humidity in the summer time of up to 50%.** Painting the edges of the coverslip is unnecessary. When the medium dries it will form a seal.

PROCEDURE

Prepare slides as required. Prior to coverslipping rinse the slides in distilled or deionized water. The excess fluid can be drained, but blotting is not recommended. Aqua-Poly/Mount is supplied in convenient 20ml dropper bottles for manual staining and in larger volumes for automated coverslippers.

Apply **two drops** of Aqua-Poly/Mount to the slide at the end or over the tissue. Carefully lower the coverslip at an angle while gently applying pressure to force any excess medium and air bubbles away from the tissue and out from under the coverslip. It is helpful to draw the coverslip to the edge of the coverslipping medium and then lower it slowly while applying pressure. Gently tilt the slide to remove any medium at the edges of the slide and coverslip. The slide can be viewed immediately or after drying. Slides can be dried at room temperature or at 4°C. Drying at 4°C will increase drying times. Do not heat the slides as this can damage or fade some stains or reactions. Slides containing fluorescent chromogens should be stored in the dark.

Aqua-Poly/Mount can be used on automated coverslippers in place of traditional solvent based coverslipping medium. The viscosity of the medium may require adjusting for proper flow rates. **Allow the liquid to come to room temperature before making any adjustments for flow rate as the viscosity will change.** The amount of medium dispensed will effect the drying time. Thicker sections will require more coverage with medium to assure the coverslip remains in place against the slide surface.

ORDERING INFORMATION

Cat #	Description	Size
18606-20	Aqua-Poly/Mount	20ml
18606-500		500ml

ADDITIONAL RELATED PRODUCTS

Cat #	Description
16864	Poly/LEM Fixative
24216	Tissue Tack Microscope Slides
22247	Poly-L-Lysine Coated Microscope Slides

TO ORDER

In The U.S. Call: 1(800) 523-2575 / (215) 343-6484
In The U.S. Fax: 1(800) 343-3291 / (215) 343-0214
In Germany Call: +(49) 6221-765767
In Germany Fax: +(49) 6221-764620
In Taiwan Call: +(49) 6221-765767
In Taiwan Fax: +(49) 6221-764620

Order online anytime: www.polysciences.com

Should any of our materials fail to perform to our specifications, we will be pleased to provide replacements or return the purchase price. We solicit your inquiries concerning all needs for life sciences work. The information given in this bulletin is to the best of our knowledge accurate, but no warranty is expressed or implied. It is the user's responsibility to determine the suitability for their own use of the products described herein, and since conditions of use are beyond our control, we disclaim all liability with respect to the use of any material supplied by us. Nothing contained herein shall be construed as a recommendation to use any product or to practice any process in violation of any law or any government regulation.