

TECHNICAL DATA SHEET 559

Page 1 of 1

BioMag® Goat anti-Rabbit IgG

Description

BioMag Goat anti-Rabbit IgG is a suspension of BioMag particles approximately 1.6µm in size, which are covalently coated with goat anti-rabbit IgG antibody. The suspension is supplied in phosphate buffered saline (pH 7.4) with EDTA and sodium azide added. After shaking vigorously or vortexing, BioMag Goat anti-Rabbit IgG is ready to use.

Characteristics

Mean Diameter: ~1.6µm
Particle Concentration: 1 mg/ml
Binding Capacity: 1mL (1mg) of BioMag Goat anti-Rabbit IgG will bind > 0.1mg of rabbit IgG.

Procedure

Researchers are advised to optimize the use of BioMag in any application as procedures designed by other manufacturers may not be ideal.

BioMag Goat anti-Rabbit IgG is suitable for use as a secondary antibody in radioimmunoassays and enzyme immunoassays that utilize a rabbit primary antibody. In general, BioMag Goat anti-Rabbit IgG may be used to quickly and conveniently separate a rabbit antibody or rabbit antibody complex from solution. (For additional information on the use of this product in cell separation, please see Technical Data Sheet 528, 'BioMag and Cell Sorting.')

Separation may be accomplished magnetically or through centrifugation.

Storage and Stability

Store at 4°C. Freezing, drying, or centrifuging BioMag may result in irreversible aggregation and loss of binding activity.

Safety

This particle suspension contains sodium azide. Sodium azide may react with lead and copper plumbing to form explosive metal azides. Upon disposal of material, flush with a large volume of water to prevent azide accumulation. Please consult the Material Safety Data Sheet for more information.

This product is for research use only and is not intended for use in humans or for *in vitro* diagnostic use.

Ordering Information

| Cat. # | Description | Size |
|-----------|------------------------------|-------|
| 84300-50 | BioMag® Goat anti-Rabbit IgG | 50ml |
| 84300-500 | BioMag® Goat anti-Rabbit IgG | 500ml |

To Order

In The U.S. Call: 1-800-523-2575 • 215-343-6484
In The U.S. FAX: 1-800-343-3291 • 215-343-0214

In Germany Call: (49) 6221-765767
In Germany FAX: (49) 6221-764620

Order online anytime at www.polysciences.com.