



U.S. Corporate Headquarters
400 Valley Road
Warrington, PA 18976
1(800) 523-2575 / (215) 343-6484
Fax: 1(800)343-3291
Email: info@polysciences.com
www.polysciences.com

Polysciences Europe GmbH
Handelsstrasse 3 D-69214
Eppelheim, Germany
+(49) 6221-765767
Fax: +(49) 6221-764620
Email: info@polysciences.de

TECHNICAL DATA SHEET 862

Page 1 of 1

IHC Hematoxylin / Bluing Kit

DESCRIPTION

IHC Hematoxylin is formulated to impart light nuclear counterstaining desired with immunohistochemical chromogen staining. Due to the controlled nuclear counterstaining the IHC Hematoxylin will not overstain and obscure the nuclear IHC staining. The counterstain and bluing solution come in one kit and are ready to use.

Volume: 500ml for A and B

PROCEDURE

1. After all IHC chromogen staining follow manufacturers recommendations for rinsing slides from any mineral oil solution and proceed with placing slides in distilled water rinse before placing in IHC ready to use hematoxylin. When counterstaining IHC slides with Hematoxylin, place in IHC hematoxylin for 1.5 minutes.
2. Rinse well in tap water 2 minutes
3. Blue in IHC Bluing Solution for 20-30 seconds
4. Rinse in tap water for 1 minute
5. Dehydrate Clear and mount

STORAGE AND HANDLING

Shipment

Storage

Store at 2° - 8° C

Safety

Refer to MSDS for details

ORDERING INFORMATION

Cat. #	Description	Size
25082-1	IHC Hematoxylin / Bluing Kit	1 kit

TO ORDER

In The U.S. Call: 1(800) 523-2575 • (215) 343-6484
In The U.S. Fax: 1(800) 343-3291 • (215) 343-0214
In Germany Call: +(49) 6221-765767
In Germany Fax: +(49) 6221-764620

Order online anytime at www.polysciences.com