



Corporate Headquarters
400 Valley Road
Warrington, PA 18976
1-800-523-2575
FAX 1-800-343-3291
Email: info@polysciences.com
www.polysciences.com

Europe - Germany
Polysciences Europe GmbH
Handelsstr. 3
D-69214 Eppelheim, Germany
(49) 6221-765767
FAX (49) 6221-764620
Email: info@polysciences.de

TECHNICAL DATA SHEET 864

Page 1 of 2

Grocott Methenamine Silver Stain (GMS) for Fungus and PCP - Room Temperature & Microwave Kit

DESCRIPTION

Grocott's alkaline hexamine-silver solution allows silver ions upon reduction, to appear dark or black, also known as "argentaaffin reaction". the ability of a silver complex solution to blacken a tissue element without the need of a reducing bath.

In fungal infections are becoming more common in the form of opportunistic infections in patients that are immunocompromised. Fungi stain weakly with hematoxylin, they are also PAS positive due to carbohydrate found in fungal cell walls.

NOTES

1. A known positive control section must be used to ensure correct differentiation has been achieved.
2. Microwave method : Volume of chromic acid must be measured
3. When microwaving solutions, ensure lids are not put on tightly to allow for expansion.
4. Prior to leaving slides to stain in microwaved solutions ensure solutions are thoroughly mixed to ensure even temperature distribution.
5. Note solutions removed from the microwave oven may spontaneously boil.
6. Times may vary according to type and power of micro-wave.
7. Combination of the original method and microwave method can be used. *ie the microwave technique for silver used with the original method for chromic acid or vice-versa.*

METHOD

1. Rehydrate sections to distilled water, use acid cleaned glassware and non metal forceps.
2. Oxidize with solution A at room temperature for 1 hour (as good time management tool, after 30 minutes, start preparing for preheated working methenamine silver solution below, allow for the waterbath to get to temperature, 60°, before making the pre-heated working solution)
3. Wash in water for a few seconds.
4. Treat sections with solution B for 1 minute
5. Wash in running tap water 3 minutes
6. Rinse thoroughly in distilled water.
7. Place in pre-heated working silver solution in a water bath at 60 °C for 15 to 20 mins until section turns yellowish-brown (Check microscopically after washing in distilled water – fungi should be dark brown).

Preparation of Preheated Working Solution

Prepare working silver solution fresh each time add in the following order in an acid cleaned coplin jar, and do not use metal forceps

Solution D ----- 23 ml
Solution C ----- 1.25 ml
Solution E ----- 3 ml
Distilled H₂O ----- 25 ml

8. Rinse thoroughly in distilled water.
9. Tone sections with solution F for 2 minutes
10. Rinse in distilled water.

8. Treat sections with solution G for 2 minutes
9. Wash with running tap water 5 minutes
10. Counterstain in working solution H 15 seconds
11. Rinse excess light green off slide with alcohol
12. Dehydrate, clear and mount

MICROWAVE METHOD

1. Bring sections to distilled water.
2. Place slides in 40 ml solution A in a loosely covered plastic coplin jar. Microwave at 150 Watt (see technical note 7) 2 min 30 sec
3. Dip slides up and down in coplin jar and leave to stand for a further 2 minutes
4. Wash in running tap water 30 seconds
5. Treat sections with solution B for 30 seconds
6. Wash in running tap water 30 seconds
7. Preheat working silver solution (loosely lidded plastic coplin jar) at 450 Watt 60 seconds

Preparation of Preheated Working Solution for Microwave

Prepare working silver solution fresh each time add in the following order in a plastic coplin jar, and do not use metal forceps, to stir solution, use an orangewood stick

Solution D ----- 23 ml
Solution C ----- 1.25 ml
Solution E ----- 3 ml
Distilled H₂O ----- 25 ml

8. Tone sections with solution F for a minute
9. Rinse in distilled water
10. Treat sections with solution G for one minute
11. Wash with running tap water 15 seconds

Working Light Green Solution Preparation

Solution H ----- 10 ml
 Distilled H₂O ----- 40 ml

12. Counterstain in working light green 15 sec
13. Rinse excess light green off slide with alcohol
14. Dehydrate, clear and mount.

RESULTS

fungi, Pneumocystis carinii, histoplasma spp ----- black
 inner parts of mycelia and hyphae ----- old rose
 leishmania spp, toxoplasma spp ----- negative
 mucin ----- dark grey
 background ----- pale green

STORAGE, SAFETY AND HANDLING

Storage

Store at room temperature

Safety

Wear protective clothing, gloves and safety glasses when preparing reagents. Refer to MSDS for more details

ORDERING INFORMATION

| Cat. # | Description | Size |
|---------|--|-------|
| 25087-1 | Grocott Methenamine Silver Stain (GMS) for Fungus and PCP - Room Temperature & Microwave Kit | 1 kit |

TO ORDER

In The U.S. Call: 1(800) 523-2575 • (215) 343-6484
In The U.S. Fax: 1(800) 343-3291 • (215) 343-0214
In Germany Call: +(49) 6221-765767
In Germany Fax: +(49) 6221-764620

Order online anytime at www.polysciences.com

Should any of our materials fail to perform to our specifications, we will be pleased to provide replacements or return the purchase price. We solicit your inquiries concerning all needs for life sciences work. The information given in this bulletin is to the best of our knowledge accurate, but no warranty is expressed or implied. It is the user's responsibility to determine the suitability for their own use of the products described herein, and since conditions of use are beyond our control, we disclaim all liability with respect to the use of any material supplied by us. Nothing contained herein shall be construed as a recommendation to use any product or to practice any process in violation of any law or any government regulation.