



U.S. Corporate Headquarters
400 Valley Rd.
Warrington, PA 18976
1(800) 523-2575 / (215) 343-6484
1(800)343-3291 fax
info@polysciences.com

Polysciences Europe GmbH
Handelsstrasse 3
D-69214 Eppelheim, Germany
+(49) 6221-765767
+(49) 6221-764620 fax
info@polysciences.de

Polysciences Asia-Pacific, Inc.
2F-1, 207 Tunhwa N. Rd.
Taipei, Taiwan 10595
(886) 2 8712 0600
(886) 2 8712 2677 fax
info@polysciences.tw

TECHNICAL DATA SHEET 848

Page 1 of 1

1X 10mM Citrate Buffer Kit, pH 6.0

for antigen retrieval FFPE, and resin

DESCRIPTION

Formalin fixation forms protein cross-links that can mask the antigenic sites in tissue specimens, thereby creating weak or false negative staining for immunohistochemical detection of certain proteins. There are currently several antigen retrieval methods commonly used and the protocols below only intended as a general guide. Polysciences, Inc. provides AR buffer kits, that include rinses and pH adjustment reagents for the 20X concentration kit. These kits are recommended for formalin fixed paraffin embedded tissues.

1. Place slides in 500ml of buffer (use a microwavable container)
2. Microwave on high for 15 minutes without a lid or cover with vented cling-film and heat at 90°C for 10 minutes (buffer must not boil)
3. Let sections stand for 15 minutes to cool
4. Wash in Tris Buffered Saline (Cat. #24082) and proceed with staining protocol

STORAGE AND HANDLING

Storage

Store at 4° C

Safety

Refer to MSDS for details

All values are considered typical based on tests believed to be accurate.
Polysciences, Inc. may change the data as appropriate.

ORDERING INFORMATION

Cat. #	Description	Size
25084	1X 10mM Citrate Buffer Kit, pH 6.0 (for antigen retrieval FFPE and resin)	1 Kit

TO ORDER

In The U.S. Call: 1(800) 523-2575 / (215) 343-6484
In The U.S. Fax: 1(800) 343-3291 / (215) 343-0214
In Germany Call: +(49) 6221-765767
In Germany Fax: +(49) 6221-764620
In Taiwan Call: +(49) 6221-765767
In Taiwan Fax: +(49) 6221-764620

Order online anytime: www.polysciences.com