


# BioMag® 15ml / 50ml Tube Separator

Catalog Number: 84102S

## DESCRIPTION

The BioMag® 15ml / 50ml Tube Separator is designed to accommodate five 15ml tubes and three 50ml tubes for a variety of applications, including cell sorting and couplings of ligands to BioMag®. Each bar magnet, embedded in a plastic rack, has a magnetic strength of 35 megagauss Oersteads.

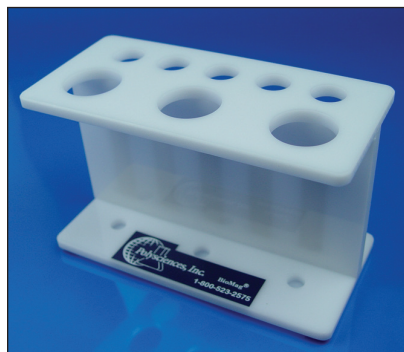

## STORAGE AND SAFETY

**Storage** The BioMag® 15ml/50ml Tube Separator should be wiped clean with mild detergent. Do not immerse. Harsh chemicals and organic solvents may damage the surface of the unit.

**Safety** Strong magnetic field. This separation unit should be used and stored a safe distance from magnetic media, such as computers, videotapes, credit cards, etc.

**This product is for research use only and is not intended for use in humans or for *in vitro* diagnostic use.**

## PROCEDURE

To use, simply place the 15ml or the 50ml tubes in the appropriate holder. The rare earth magnets will draw the BioMag® solution toward the wall of the separator to complete separation in 3 to 10 minutes. For 15ml tubes, separations are optimized when volumes of BioMag® solution are between 7ml and 12ml. For 50ml tubes, separations are optimized when volumes of BioMag® solution are between 25ml and 40ml. Once a separation is complete, aspirate the supernatant without disturbing the separated BioMag®.

## ORDERING INFORMATION

Cat. #	Description	Size
84102S-1	BioMag® 15ml / 50ml Tube Separator	1 each

## TO ORDER

In The U.S. Call: 1(800) 523-2575 • (215) 343-6484  
In The U.S. Fax: 1(800) 343-3291 • (215) 343-0214

In Germany Call: +(49) 06201-845200  
In Germany Fax: +(49) 06201-8452020

In Asia Call: (886) 2 8712 0600  
In Asia Fax: (886) 2 8712 2677

Order online anytime at [www.polysciences.com](http://www.polysciences.com)