



U.S. Corporate Headquarters
400 Valley Rd.
Warrington, PA 18976
1(800) 523-2575 / (215) 343-6484
1(800)343-3291 fax
info@polysciences.com

Polysciences Europe GmbH
Badener Str. 13
69493 Hirschberg an der Bergstrasse,
Germany
+(49) 06201-845200
+(49) 06201-8452020 fax
info@polysciences.de

Polysciences Asia-Pacific, Inc.
2F-1, 207 DunHua N. Rd.
Taipei, Taiwan 10595
(886) 2 8712 0600
(886) 2 8712 2677 fax
info@polysciences.tw

TECHNICAL DATA SHEET 574

Page 1 of 1

BioMag® Multi-32 Microcentrifuge Tube Separator

Catalog Number: 84106S

DESCRIPTION

The BioMag® Multi-32 Microcentrifuge Tube Separator holds thirty-two microcentrifuge tubes for molecular biology and other applications. Sixteen tubes may be magnetically separated at a time.

PROCEDURE

Magnetic separations can be performed on up to sixteen microcentrifuge tubes simply by placing the tubes in the inner row of the separator. Sixteen tubes may also be held in the outer rows of the separator for mixing and short incubations. Holds 1.0ml and 1.5ml microcentrifuge tubes. The BioMag® solution will move toward the side of the separator in less than 5 minutes to complete the separation. Pellets will form on the side of the test tube. Once separation is complete, aspirate the supernatant with the BioMag® Separator in place. Be careful not to disturb the separated BioMag®.

STORAGE AND SAFETY

Storage The BioMag® Multi-32 Microcentrifuge Tube Separator should be wiped clean with mild detergent. Do not immerse. Harsh chemicals and organic solvents may damage the surface of the unit.

Safety Strong magnetic field. This separation unit should be used and stored a safe distance from magnetic media, such as computers, videotapes, credit cards, etc.

This product is for research use only and is not intended for use in humans or for *in vitro* diagnostic use.



ORDERING INFORMATION

Cat. #	Description	Size
84106S-1	BioMag® Multi-32 Microcentrifuge Tube Separator	1 each

TO ORDER

In The U.S. Call: 1(800) 523-2575 • (215) 343-6484

In The U.S. Fax: 1(800) 343-3291 • (215) 343-0214

In Germany Call: +(49) 06201-845200

In Germany Fax: +(49) 06201-8452020

In Asia Call: (886) 2 8712 0600

In Asia Fax: (886) 2 8712 2677

Order online anytime at www.polysciences.com

Should any of our materials fail to perform to our specifications, we will be pleased to provide replacements or return the purchase price. We solicit your inquiries concerning all needs for life sciences work. The information given in this bulletin is to the best of our knowledge accurate, but no warranty is expressed or implied. It is the user's responsibility to determine the suitability for their own use of the products described herein, and since conditions of use are beyond our control, we disclaim all liability with respect to the use of any material supplied by us. Nothing contained herein shall be construed as a recommendation to use any product or to practice any process in violation of any law or any government regulation.