

**U.S. Corporate Headquarters**  
400 Valley Rd.  
Warrington, PA 18976  
1(800) 523-2575 / (215) 343-6484  
1(800)343-3291 fax  
info@polysciences.com

**Polysciences Europe GmbH**  
Handelsstrasse 3  
D-69214 Eppelheim, Germany  
+(49) 6221-765767  
+(49) 6221-764620 fax  
info@polysciences.de

**Polysciences Asia-Pacific, Inc.**  
2F-1, 207 Dunhua N. Rd.  
Taipei, Taiwan 10595  
(886) 2 8712 0600  
(886) 2 8712 2677 fax  
info@polysciences.tw

## TECHNICAL DATA SHEET 119A

Page 1 of 2

# Osmium Tetroxide, 99.95%

### PACKAGING:

All of our osmium tetroxide is packaged in pre-scored unlabeled ampoules. Each ampoule of solution is packaged with an ampoule cracker.

### BACKGROUND AND PROPERTIES:

Although the material dissolves slowly in water; ~7.24% at 25°C solubility, it is 518 times more soluble in CCl<sub>4</sub> (carbon tetrachloride).

- Volatilizes readily at room temperature.
- Vapor pressure is given by  $\log p = 10,000/4.57T + 5.49$

A 1% solution of OsO<sub>4</sub> in distilled water is 0.04 mol/L.<sup>1</sup> Aqueous solutions of OsO<sub>4</sub> are hypotonic. To increase osmolarity of the final fixing solution, electrolytes or non-electrolytes are used.<sup>1</sup>

### PREPARATION OF OsO<sub>4</sub> SOLUTIONS:

Polysciences' ampoules of OsO<sub>4</sub> are offered in both crystalline and solution form. For those wishing to have aqueous solutions available in fresh uncontaminated form, we suggest our vials containing 4% aqueous OsO<sub>4</sub>. This can be reduced to 2% or 1% OsO<sub>4</sub> by adding the 4% solution to an appropriate volume of the desired buffer. A more economic form is our high purity crystalline OsO<sub>4</sub> which can be easily dissolved.

A stock solution of Osmium for buffering is prepared from crystalline as follows: The sealed ampoule and the storage bottle are thoroughly cleaned and rinsed. Gloves should be worn at this step not only for handling purposes, but also to prevent any organic contamination of the ampoule by direct skin contact.

The 2% OsO<sub>4</sub> is made by opening the ampoule containing 1.0 gram OsO<sub>4</sub> and placing the ampoule into 50 ml of glass-distilled water in a clean storage vessel. This solution is stable and keeps at 4°C for many months if kept clean and tightly closed. OsO<sub>4</sub> crystals will dissolve more rapidly using a magnetic stirrer, shaker, or sonicator. Suitable buffer systems can be mixed in any proportion with a stock solution of osmium tetroxide to give the desired concentration of fixative and buffer.

Osmium tetroxide in carbon tetrachloride solution has been used to fix algae. The fixative diffuses from the non-aqueous to the aqueous phase and provides a condition similar to vapor fixation but at a higher concentration of OsO<sub>4</sub>.<sup>2</sup>


### PRECAUTIONS AND FIRST AID:

OsO<sub>4</sub> vapors and solutions stain skin black and are very toxic and damaging to the eyes, skin and respiratory tract. Do not breathe vapors. The effects of contact may be delayed. This material is a strong oxidizer. Wear goggles and gloves when working with OsO<sub>4</sub>. Handle material under a hood.

**Eyes:** Flush with plenty of water for at least 15 minutes. Get prompt medical attention.

**Skin:** Wash with diluted sodium bisulfite solution, followed by soap and water.

**Inhalation:** Remove patient to fresh air; if not breathing, give artificial respiration. Get prompt medical attention. Remove contaminated clothing.

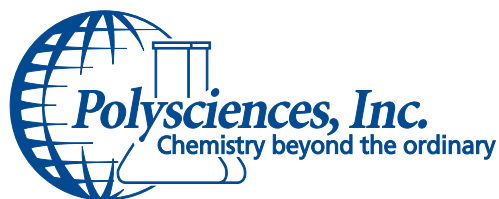

**U.S. Corporate Headquarters**  
400 Valley Rd.  
Warrington, PA 18976  
1(800) 523-2575 / (215) 343-6484  
1(800)343-3291 fax  
info@polysciences.com

**Polysciences Europe GmbH**  
Handelsstrasse 3  
D-69214 Eppelheim, Germany  
+(49) 6221-765767  
+(49) 6221-764620 fax  
info@polysciences.de

**Polysciences Asia-Pacific, Inc.**  
2F-1, 207 Dunhua N. Rd.  
Taipei, Taiwan 10595  
(886) 2 8712 0600  
(886) 2 8712 2677 fax  
info@polysciences.tw

## TECHNICAL DATA SHEET 119A

Page 2 of 2

### ORDERING INFORMATION:

Cat. #	Description	Size
0223A	Osmium tetroxide, crystalline 99%	5 x 1g
0223B	Osmium tetroxide, crystalline 99%	10 x 1g
0223C	Osmium tetroxide, crystalline 99%	10 x 1/2g
0223D	Osmium tetroxide, crystalline 99%	10 x 1/4g
0972A	Osmium tetroxide, 4% microfiltered solution	20 x 2mL amp
0972B	Osmium tetroxide, 4% microfiltered solution	5 x 10mL amp
0972C	Osmium tetroxide, 4% microfiltered solution	20 x 10mL amp
23310	Osmium tetroxide, 2% microfiltered solution	10 x 2mL amp
23311	Osmium tetroxide, 2% microfiltered solution	10 x 5mL amp

### REFERENCES:

1. Hayat, M.A., Fixation for Electron Microscopy, Academic Press, 148 (1981), available from Polysciences, Inc. (Cat. No. 17459)
2. Hobbs, M.J., Stain Tech., **44**, 217 (1969).

### TO ORDER

In The U.S. Call: 1(800) 523-2575 • (215) 343-6484  
In The U.S. Fax: 1(800) 343-3291 • (215) 343-0214

In Germany Call: +(49) 6221-765767  
In Germany Fax: +(49) 6221-764620

In Asia Call: (886) 2 8712 0600  
In Asia Fax: (886) 2 8712 2677

Order online anytime at [www.polysciences.com](http://www.polysciences.com)